Thermo Scientific LipidSearch Software for Lipidomics Workflows

Automated Identification and Relative Quantitation of Lipids by LC/MS





The promise of lipidomics

Lipidomics is a new field of study crucial for understanding cellular physiology and pathology. The application of lipidomic profiling to disease phenotype analysis is a rapidly growing aspect of translational medical research. Identification of unique lipid biomarkers has the potential to distinguish healthy individuals from individuals at risk for disease, detect diseases earlier, and facilitate development of personalized treatments.

Liquid chromatography combined with mass spectrometry (LC/MS) is a widely adopted technique for lipidomics analyses. Relative and absolute quantitation, and identification, of lipids from biological samples requires sophisticated software with an extensive, comprehensive database. Thermo Scientific[™] LipidSearch[™] software provides accurate identification of lipids and automatically integrates complex data into a concise report. With its easy-to-use web-based interface, it dramatically reduces data analysis time.

A	utomated	identi	fic
q	uantitatio	n with	Lip

LipidSearch software, developed jointly by Professor Ryo Taguchi and MKI, (Tokyo, Japan), is a powerful new tool for automatic identification and relative quantification of cellular lipid molecular species from large amounts of mass spectrometric data obtained in nano-infusion or LC-MS experiments. Using the industry-leading, high-resolution, accurate-mass Thermo Scientific Orbitrap[™] technology with exclusive LipidSearch software, the most accurate and confident lipid profiles and identifications can be achieved more quickly than ever before.

- Compatible with data acquired from Thermo Scientific[™] triple quadrupole, ion trap, and Orbitrap mass spectrometers
- Largest lipid database containing >1.5 million lipid ions and their predicted fragment ions
- Identification algorithms for product ion, precursor ion and neutral loss scans
- Alignment of lipid data obtained from multiple LC-MS and MSⁿ experiments
- Relative quantitation of identified lipid precursors in either LC-MS or infusion experiments

The database contains defined structures and includes more than **1.5 million lipid ions and their** predicted fragment ions. Fragmentation patterns are calculated and improved by using expert knowledge based on experimental results. Lipid adduct ions and MSⁿ fingerprints are also included. Data are stored in XML files and are easily customized.

F	atty acid	ls		
sn1 CH₂	O II —OC-R ₁			
			Adduct io	n
sn2 CH ₂ —	O II —OC-R ₂		о НС-0 ⁻	N
sn3 CH₂	0 	(-O-CH ₂ -CH ₂ N - - (CH₃ I≐CH₃ CH₃	
Lipid backbone	Lip	oid headgr	oup	- F

nit a Explore									
	r P	arameters							
c:\lipidsearch\	data -	Identification	Quantitation	Filters	Class	Ion			
data	1.000	Batch			Search op	tions			
	st_1.raw st_1.raw	Job Name	YEAST LIPIDS		Save o	ptions	ProductSearch_	QEX	~
AKOPost	_1.raw _1.raw	Comment	1		Search Exp1	туре Туре	Parent LC-MS	Product O 7	arget 🕘 Mult
A2-top5-	neg.raw),raw Database		Parer	t tol	0.1	Da 💌		
B2-top5-	neg.raw			NL/Pr	ec tol	0.5	Da 🔹		
C2-top5-	5-neg.raw 5-neg.raw Target Databa _Neg.raw	General C Exactive		Precur	sor tol	5.0	ppm -		
StdMix_M		Orbitrap		Produ	ct tol	8.0	ppm •		
StdMix_F	Pos.raw		0.0.012		Merge Rar	nge(min.)	99999.0		
clear		Peak detect	ion		Min Peak W	/idth(min.)	0.0		
		Recalc Isote	ope 💿 on 😳 off				parent 0.01		
		R.T. interval(Min) 0.0		Intensity	threshold	threshold type		
		R.T. Range(I	Hin)				product ion 5.0	absolute (%)	
					m-Score I	breshold	3.0		

Easy-to-use web browser interface for automated and reliable identification

ation and relative pidSearch software



Lipid fragment ions related to lipid headgroup, fatty acids, and backbone





From peak detection to relative quantitation and identification, LipidSearch software provides an easy-to-use, automated workflow.



Smart peak detection

Step 2

Identification

Scoring algorithms

• Group-specific ID (targeted)

• Comprehensive ID (untargeted)

The peak detection engine implemented in LipidSearch software can handle different MS experiments and platforms. The combination of unique peak detection algorithms, tailored for each experiment and instrument type, and mass spectral processing functions enables accurate peak detection.

loss scans from lipids mixtures. Soli Name Test Identification Module - Lipid

lipids based on the polar head groups or fatty acids using a combination of

precursor ion scanning and neutral

stored in the database.

• LipidSearch software also provides a set of scoring algorithms to filter out lower probability results.



C. Relative quantitation by lipid class



Accurate peak areas are computed by denoising and smoothing the peak profiles prior to separating any partially overlapped peaks. Comparative analysis is then carried out between the multiple sample and control groups using *t*-test statistics. The mean peak area result for each group is displayed in a "box and whisker" plot.



Alignment results for CoQ_7 , $[M+NH_4]^+$



Lipid profiling of wild type and knockout yeast strains reveals detailed changes for individual lipid species

Wild-type (WT) yeast (S. Cerevisiae) continue to grow after glucose is exhausted from the media (diauxic shift point) whereas knockout (KO) yeast have a defect in CoQ production and do not grow after the shift point. Duplicates of WT and KO yeast were collected post shift for lipidomics LC-MS analysis using a Thermo Scientific[™] Q Exactive[™] hybrid quadrupole-Orbitrap MS both in MS and data-dependent MS² mode.

(Peake, D.A.; Wang, J.; Huang, P.; Jochem, A.; Higbee, A.; Pagliarini, D.J., Quantitative yeast lipidomics via LC-MS profiling using the Q Exactive Orbitrap mass spectrometer, presented at the LIPID MAPS Annual Meeting 2012, May 7-8, 2012, La Jolla, CA.)





Data Processing

LC-MS raw files containing full scan MS and data-dependent - MS² data were searched for FA (fatty acids), sphinganine, Lyso-GP (glycerophospholipids), MG (mono-acyl glycerol), GP (PA, PC, PE, PG, PI, PS), Cer (ceramides) and CoQ (co-enzyme) lipid classes using a mass tolerance of 5 ppm for precursor ions and 8 ppm for product ions. The search results from 2 WT and 2 KO samples were aligned using a 0.25 min tolerance window and a combined report was generated. A total of 738 lipid isomers with 542 different formulas were identified and correlated between the 4 data files.



LC-MS chromatogram of lipids from yeast

	Name	Eļ.	RawData	🗐, Туре	🗐 ЕхрТуре	Proce	55	Result
Yeast ①	1	11WTPost_1.raw		Product	LC	PIQ	٢	8322 2639
Yeast ()	3	12WTPost_1.raw		Product	LC	PIQ	٢	8332 2638
Yeast ()	3	3KOPost_1.raw		Product	LC	P I Q	٢	8309 2637
Yeast ①)	4KOPost_1.raw		Product	LC	PIQ	•	8306 26366

Aligned Results Summary

E. Na	me	🗐 Туре	🗐 ЕхрТуре	EI, P	rocess	🗐 Result
Yeast (All)	0	Product	LC	м	٢	88738 2814
Yeast (Ma	in) ①	Product	LC	М	(1)	8 542 8 607

LipidSearch status view of search and alignment results

Step 2

Identification

For each MS² spectrum, search results are summarized, along with a score indicating the fit, for lipid species matching the predicted fragmentation pattern from the database. If a mixture of lipids is found, the most abundant lipid is displayed. The fragment ions used to identify the lipid species are highlighted in red when the match is selected.

ID of PE 16:0-18:1, MS² of m/z 716.5230 [C₃₉H₇₅NO₈P]⁻



Step 3

Alignment and Quantitation

Alignment results are summarized for annotated lipid species eluting within the retention time tolerance window. The relative peak areas are reported along with the *t*-test statistics. The aligned chromatograms for two Lyso - PC 18:1 isomers are shown to the right. The structural assignments are based on the known elution order of lyso-phospholipids on reversed-phase HPLC columns















Using the industry-leading, high-resolution, accurate-mass Thermo Scientific[™] Orbitrap[™] technology with exclusive LipidSearch software, achieve the most accurate and confident lipid profiles and identifications more quickly than ever before. For targeted quantitation, LipidSearch software also supports the Thermo Scientific TSQ triple quadrupole MS systems. For both lipidomics workflows, LipidSearch software automatically integrates complex data into a combined report to dramatically reduce data analysis.

Minimum/Recommended PC Requirements

OS: Microsoft® Windows® 7 Professional (x64) or Windows 8 (x64); English language
CPU: Quad - or multi-core CPU, 2 GHz or higher
Memory: 8 GB RAM or higher
HDD: 50 GB free space for LipidSearch software; 500 GB hard drive
SSD (optional): 50 GB free space for LipidSearch software; 256 GB solid-state drive



For Research Use Only. Not for use in diagnostic procedures.

www.thermoscientific.com

©2014 Thermo Fisher Scientific Inc. LipidSearch is a registered trademark of Mitsui Knowledge Industry Co., LTD. Microsoft and Windows are registered trademarks of Microsoft Corp. All rights reserved. ISO is a trademark of the International Standards Organization. All other trademarks are the property of Thermo Fisher Scientific and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.



Singapore +65 6289 1190 Spain +34 914 845 965 Sweden +46 8 556 468 00 Switzerland +41 61 716 77 00 UK +44 1442 233555 USA +1 800 532 4752



Africa +43 1 333 50 34 0 Australia +61 3 9757 4300 Austria +43 810 282 206 Belgium +32 53 73 42 41 Canada +1 800 530 8447 China 800 810 5118 (ree call domestic) 400 650 5118 BR63951 F 01/14

Denmark +45 70 23 62 60 Europe-Other +43 1 333 50 34 0 Finland +358 9 3291 0200 France +33 1 60 92 48 00 Germany +49 6103 408 1014 India +91 22 6742 9494 Italy +39 02 950 591 Japan +81 45 453 9100 Latin America +1 561 688 8700 Middle East +43 1 333 50 34 0 Netherlands +31 76 579 55 55 New Zealand +64 9 980 6700 Norway +46 8 556 468 00 Russia/CIS +43 1 333 50 34 0