

Mass spectrometry

TSQ Altis Plus EFOX Edition Triple Quadrupole Mass Detector

Complex contaminant analysis redefined

Benefits

- Dual-CID expands traditional MS² method development for greater flexibility, stronger confidence, and simplified optimization
- QR5 Plus segmented quadrupoles with hyperbolic surfaces allow running high-resolution selective-reaction monitoring (H-SRM) transitions up to 0.2 Da FWHM for even more consistency, accuracy, reliability, and reproducibility
- Industry-leading selected-reaction monitoring (SRM) speeds enable quantitation of more compounds in less time
- Provides robustness, reliability, and consistency that are unprecedented in a high-end triple quadrupole mass spectrometer
- Gives industry-leading polarity switching times that include signal stabilization
- Provides automated compound optimization and intuitive instrument interface
- Access to the comprehensive Thermo Scientific™ mzCloud™ compound database

The Thermo Scientific™ TSQ Altis™ Plus EFOX Edition Mass Detector redefines ultimate performance and flexibility to address demanding quantitation workflows with ease. With superior acquisition speeds, enhanced sensitivity, and stability, the TSQ Altis Plus EFOX Edition MS delivers unprecedented accuracy and precision for low-level compounds in complex matrices. Advanced control software, database integration, and templates streamline rigorous method development. Combined with powerful Thermo Scientific™ chromatographic systems, and leading, fit-for-purpose Thermo Scientific™ Chromeleon™ Chromatography Data System (CDS) software, a TSQ Altis Plus EFOX Edition solution provides the highest performance for your most challenging quantitative applications.




Hardware features

Thermo Scientific™ OptaMax™ NG API source

- Automatic connection of all gases and voltages on installation simplifies operation and improves reliability
- Automatic source recognition simplifies use and data logging
- Sweep gas reduces chemical noise
- Enhanced exhaust port efficiently removes solvent vapor, improving uptime and reducing chemical noise
- Flexible X, Y, and Z positioning for all ionization probes maximizes performance
- Optimal 60-degree spray angle

TSQ Altis Plus EFOX Edition ion flight path

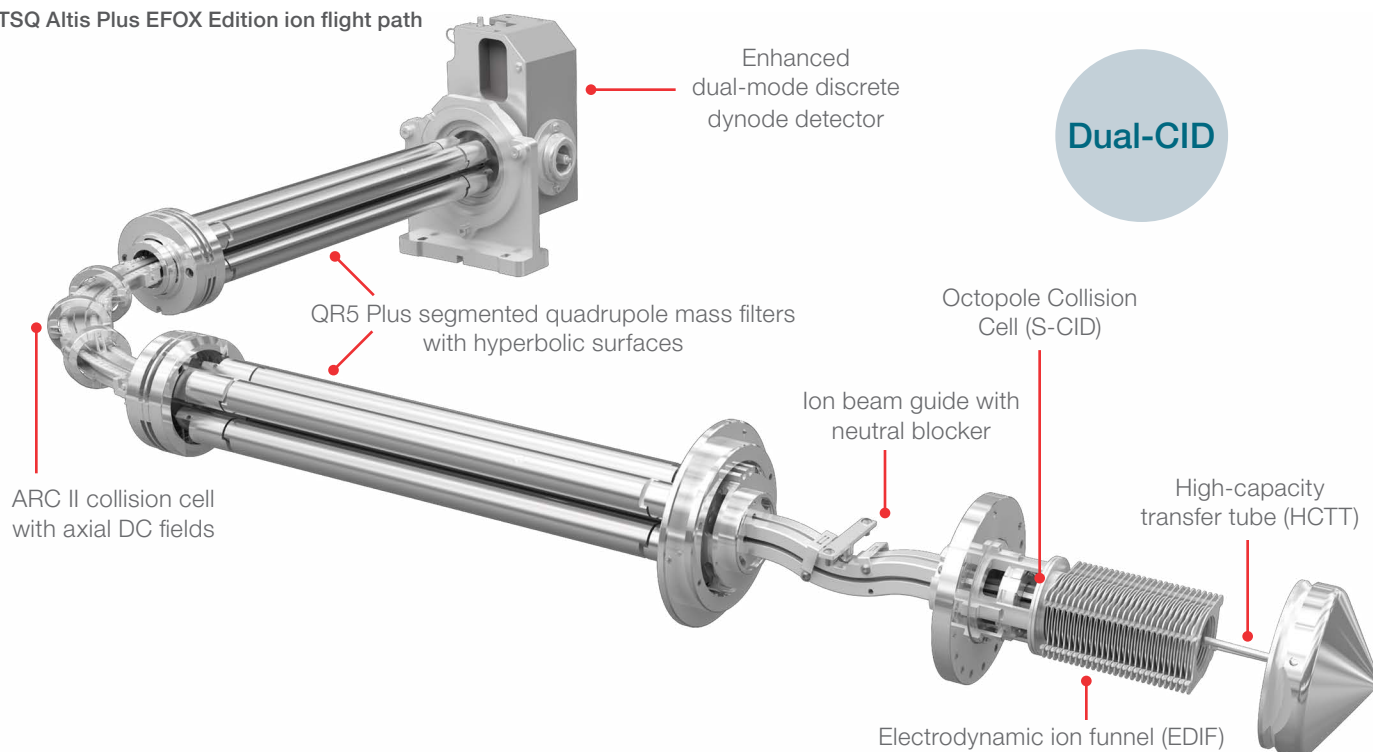

- Integrated APCI functionality with interchangeable HESI and APCI ionization probes
- APPI compatibility
- Dedicated mass calibration probe

High-capacity transfer tube

The optimized high-capacity transfer tube (HCTT) transfers more ions into the vacuum system for improved sensitivity while maintaining robustness and ease of use.

Ion optics

Electrodynamic ion funnel RF lens

The electrodynamic ion funnel (EDIF) efficiently captures ions as they leave the transfer tube. Its broad transmission curve reduces ion losses, increasing sensitivity. By design, the EDIF gently moves ions from atmosphere to vacuum, reducing in-source fragmentation.

Octopole Collision cell (S-CID)

The S-CID enables pre-quadrupole CID fragmentation using nitrogen. It offers a similar fragmentation pattern as ARC II cell without any parent ion selection. The combination of the S-CID and ARC-CID (Dual-CID fragmentation) allows generation of MS³-like fragment ions.

Ion beam guide and neutral blocker

The ion beam guide, with its neutral blocker, stops neutrals and high-velocity clusters, keeping the ion path cleaner, reducing noise, and increasing sensitivity and robustness.

QR5 Plus segmented quadrupole mass filters with hyperbolic surfaces (Q1 and Q3)

Segmented hyperbolic-surface quadrupole mass filters with 5.25 mm field radius deliver industry-leading ion transmission across the mass range at resolutions up to 0.2 Da FWHM with improved mass stability for the ultimate in sensitivity at all masses and resolutions.

Active Reaction Collision cell II (ARC II)

The 90-degree, high-pressure argon-filled collision cell produces efficient fragmentation for high sensitivity, while guiding ions away from neutral reaction products to reduce noise. The new design improves stability and transmission of low-mass product ions as well as product ions generated from large m/z differences between precursor and product ions, critical for key applications. The axial DC field speeds ion transits through the collision cell yielding up to 600 SRMs/sec with zero cross-talk, providing excellent sensitivity and speed.

Detector

- Discrete-dynode detector with increased surface area greatly extends the detector lifetime
- Dual-mode function increases sensitivity by operating in pulse-counting mode when ion flux is low and analog mode when ion flux is high
- Greater than six orders of dynamic range provide high-confidence quantitation

Vacuum system

- Four-stage differentially pumped vacuum manifold
- Advanced triple-inlet turbomolecular pump integrated with the vacuum manifold
- Environmentally friendly dry-pump option available

Integrated syringe pump

Fully automated data system control of the syringe pump with plug-and-play support.

Optional divert valve

Automated control of the divert valve enables switching of the solvent front, gradient end point, or any portion of the HPLC run to waste.

Optional ion sources

- Thermo Scientific™ OptaMax™ Duet NG API ion source enables HESI and APCI functionality without manual changes in the hardware configuration

Software features

Data system

- High-performance PC with Intel™ microprocessor
- High-resolution LCD color monitor
- Microsoft™ Windows™ 10 operating system (64-bit)

Standard MS software

- Chromeleon CDS software unifies workflows for routine food and environmental quantitative mass spectrometry analysis
- Tune editor for system calibration, diagnostics, and manual data acquisition
- Fully automated loop injection or infusion-based compound optimization directly incorporated into the instrument method
- Method editor with comprehensive application-specific template library and intuitive user interface to facilitate method development
- Thermo Scientific™ mzCloud™ database access to allow importation of compound name, SRM transitions, and collision energies
- Dwell time weighting factor for optimizing efficient SRM acquisition
- Visual display of relationship between chromatographic peak width and data points per peak to determine optimum quantitative performance

Scan functions

- Highly sensitive full-scan mass spectrometry
- Selected-ion monitoring (SIM)
- SRM with up to 30,000 SRMs definable and up to 800 SRMs/sec and timed SRMs

- Dual Stage Reaction Monitoring (DSRM) and timed DSRMs
- High-resolution (0.2 Da FWHM) selected-reaction monitoring (H-SRM)
- Product ion scan
- Precursor ion scan
- Neutral-loss scan
- Reverse energy ramp MS/MS spectra (RER) gives information-rich MS/MS spectra for compound identification
- Polarity switching capabilities
- Quantitation-enhanced data-dependent (QED) MS/MS
- Mixed Mode Scan Function
- Automated on-the-fly adjustment of retention time windows to accommodate chromatographic fluctuations
- mzCloud database integration allows access to over 19,000 compounds

Performance specifications

Sensitivity*

Positive electrospray (HESI)

A 5 µL injection of a 200 fg/µL reserpine solution will produce a minimum signal-to-noise ratio of 1,500,000:1 for the transition of the protonated molecule at m/z 609.3 to the fragment ion at m/z 195.1 when operated in SRM mode with Q1 and Q3 resolution set to 0.2 and 0.7 Da FWHM respectively. The IDL for the TSQ Altis Plus EFOX Edition MS is <0.2 fg, calculated from 10 injections of 0.5 fg reserpine on the column.

Atmospheric pressure chemical ionization (APCI)

A 5 µL loop injection of a 200 fg/µL reserpine solution will produce a minimum signal-to-noise ratio of 100,000:1 for the transition of the protonated molecular ion at m/z 609.3 to the fragment ion at m/z 195.1 when operated in SRM mode with Q1 and Q3 resolution both set to 0.7 Da FWHM.

Negative electrospray (nHESI)

A 5 µL loop injection of a 200 fg/µL chloramphenicol solution will produce a minimum signal-to-noise ratio of 1,500,000:1 for the transition of the deprotonated molecular ion at m/z 321.0 to the fragment ion at m/z 152.0 when operated in SRM mode with Q1 and Q3 resolution set to 0.7 Da FWHM. The IDL for the TSQ Altis Plus EFOX Edition MS is <0.25 fg, calculated from 10 injections of 0.6 fg chloramphenicol on the column.

Mass range

m/z 2–2010

Resolution

Q1 and Q3 adjustable to 0.2 Da peak width (FWHM) across the entire mass range

Scan rate

- 15,000 amu/sec at a resolution of 2 Da FWHM
- Up to 800 SRMs/sec (for resolutions from 0.2 through 2.0 FWHM)
- Polarity switching: 5 msec electronic switching time that includes signal stabilization

Mass stability

Mass assignment will be within ± 0.1 Da over a 24-hour period. The laboratory room temperature must be maintained between 18–27 °C (65–81 °F). The room temperature may not change by more than 5 °C (9 °F) during this period.

Installation requirements

Power

- Three 230 Vac $\pm 10\%$, 50/60 Hz at 16 A minimum
- Four 120 Vac ± 6 –10%, 50/60 Hz at 20 A or four 230 Vac $\pm 10\%$, 50/60 Hz at 13 A

Gas

- Collision gas: 99.995% pure Argon (99.995% Nitrogen also supported)
- Collision gas supply pressure: 135 ± 70 kPa (20 ± 10 psig)
- Sheath/aux/sweep gas: 99% pure Nitrogen
- Sheath/aux/sweep gas supply pressure: 690 ± 140 kPa (100 ± 20 psig)
- Maximum gas consumption: ~20 L/min
- Maximum gas consumption with the Thermo Scientific™ FAIMS Pro Duo interface installed: ~55 L/min

Environment

- Functional temperature range: 15–27 °C (59–81 °F)
- Optimal temperature range: 18–21 °C (65–70 °F)
- MS Power consumption in scan mode (Heat output): 1,100 W (3,740 Btu/h)
- Total system power consumption (Heat output) with dry pump: 1,690 W (5750 Btu/h)
- Total audible noise for TSQ Altis Plus EFOX Edition MS with dry pump: 55 dB

Dimensions

Size

- TSQ Altis Plus EFOX Edition MS: 680 × 760 × 840 mm (H × W × D – 27 × 30 × 33 in.)
- Edwards™ nXR120i dry forepump: 301 × 217 × 494 mm (H × W × D – 11.8 × 8.5 × 19.4 in.)

Weight

- TSQ Altis Plus EFOX Edition MS: 125 kg (275 lb)
- Edwards™ nXR120i dry forepump: 29 kg (64 lb)

*The signal-to-noise ratio is a manufacturing specification only. The signal-to-noise ratio specifies the minimum system response expected in accordance to the performance test for reserpine and chloramphenicol evaluated under the Thermo Fisher Scientific guidelines. The specified signal-to-noise ratio in the product specification sheet does not represent the limit of detection or limit of quantitation of the mass spectrometer or any assay. The TSQ Altis Plus EFOX Edition MS is tested and installed according to the TSQ Altis Plus EFOX Edition Installation Work Instructions using the Small Molecule Standard Solution. With the new installation procedure, we will officially withdraw from running the Signal-to-Noise sensitivity specs during the system installation. Signal to noise data collected at our manufacturing facility will be available to review upon request. TSQ Altis Plus system is tested and installed according to TSQ Altis Plus MS installation Work Instructions.

 Learn more at thermofisher.com/efox